

HISTOLOGY PEN



| CODE | DESCRIPTION | Packaging |
|-------|------------------------------|-----------|
| 11-50 | Pen for cassettes and slides | 12 pcs |

Distributor: Bio-Optica Milano S.p.A.

Date of issue: 06/09/2018
Rev. 01

TECHNICAL FEATURES

Permanent marking pen for writing on plastic cassettes and glass slides.

Ideal in histology and cytology labs to classify cassettes and slides.

The ink resists to the chemical agents more used in the labs, like for example formalin, alcohols and xylene (and all its substitutive).

The writings remain clean and legible also after all labs procedures, including processing and staining procedures. The fine tip allows writing clearly.

The ink dries fast, without leaving any dribble.

Ink colour: black.

Point thickness: about 1 mm.

Writing tip material: rigid nylon

Mode of use: uncap the pen and start writing. The ink will be dried in about 30 seconds.

To avoid running, do not touch the ink with wet hands or dip in water before it is dried.

Store in vertical position, as indicated on the packaging.

If correctly closed, the product has a validity of 5 years.

Product for generic laboratory uses only. It does not request a CE mark.

Date of issue: 06/09/2018

Rev. 01